Instruction Manual Mini Auto Stirrer

EN - English		1
FR - Français		7
ES - Español		14
IT - Italiano		21
DE - Deutsch		28
PT - Português		35
NL - Nederlands		42
NO - Norsk		46
DA - Dansk		50
SV - Svenska		54
FI - Suomi		58
HU - Magyar		62
PL - Polski		66

European Catalogue Numbers:

Euro Plug: 444-2856

UK Plug: 444-2857

Swiss Plug: 444-2858

TABLE OF CONTENTS

Package Contents	1
Warranty	1
Installation	2
Maintenance & Servicing	2
Intended Use	2
Equipment Disposal	2
Environmental Conditions	2
Safety Instructions	3
Warning	3
Standard & Regulations	3
Specifications	4
Operating Instructions	5
Technical Service	6
Troubleshooting	6
EC Declaration of Conformity	70-71

PACKAGE CONTENTS

Mini Auto Stirrer
 Stir bar
 183cm power cord
 Instruction manual

WARRANTY

VWR International warrants that this product will be free from defects in material and workmanship for a period of two (2) years from date of purchase. If a defect is present, VWR will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear. For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products. IT IS EXPRESSLY AGREED THAT THIS WARRANTY WILL BE IN LIEU OF ALL WARRANTIES OF FITNESS AND IN LIEU OF THE WARRANTY OF MERCHANTABILITY.

Legal Manufacturer:

VWR International bvba • Researchpark Haasrode 2020 • Geldenaakesbaan 464 • B-3001 Leuven • + 32 16 385011 • <http://be.vwr.com>

INSTALLATION

Upon receiving the VWR Mini Auto Stirrer, check to ensure that no damage has occurred during shipment. It is important that any damage that occurred in transport is detected at the time of unpacking. If you do find such damage the carrier must be notified immediately.

After unpacking, place the Mini Auto Stirrer on a level bench or table, away from explosive vapors. Ensure that the surface on which the unit is placed will withstand typical heat produced by the unit. Always place the unit on a sturdy work surface.

The Mini Auto Stirrer is supplied with a 3-prong power cord that should be plugged into a standard 3-prong grounded outlet. If the cord supplied does not meet your needs please use an approved power cord that suits local codes and electric supply. Replacement of the plug must be made by a qualified electrician.

MAINTENANCE & SERVICING

The Mini Auto Stirrer is built for long, trouble-free, dependable service. No lubrication or other technical user maintenance is required. The unit should be given the care normally required for any electrical appliance. Avoid wetting or unnecessary exposure to fumes. Spills should be removed promptly.

DO NOT use a cleaning agent or solvent on the front panel which is abrasive or harmful to plastics, nor one which is flammable. Always ensure the power is disconnected from the unit prior to any cleaning. If the unit ever requires service, contact your VWR representative.

INTENDED USE

The Mini Auto Stirrer is intended for general laboratory use.

EQUIPMENT DISPOSAL

This equipment must not be disposed of with unsorted waste. It is your responsibility to correctly dispose of the equipment at life-cycle-end by handing it over to an authorized facility for separate collection and recycling. It is also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect the persons involved in the disposal and recycling of the equipment from health hazards.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment. By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

ENVIRONMENTAL CONDITIONS

Operating Conditions: Indoor use only.

Temperature:	5 to 40°C
Humidity:	20% to 85% relative humidity, non-condensing
Altitude:	0 to 2000 M above sea level

Non-Operating Storage:

Temperature:	-20 to 65°C
Humidity:	20% to 85% relative humidity, non-condensing

Installation Category II and Pollution Degree 2 in accordance with IEC 664.

SAFETY INSTRUCTIONS

Please read the entire instruction manual before operating the Mini Auto Stirrer.

WARNING! DO NOT use the Mini Auto Stirrer in a hazardous atmosphere or with hazardous materials for which the unit was not designed. Also, the user should be aware that the protection provided by the equipment may be impaired if used with accessories not provided or recommended by the manufacturer, or used in a manner not specified by the manufacturer.

Always operate unit on a level surface for best performance and maximum safety.

DO NOT lift unit by the top plate.

CAUTION! To avoid electrical shock, completely cut off power to the unit by disconnecting the power cord from the wall outlet. Disconnect unit from the power supply prior to maintenance and servicing.

Spills should be removed promptly. **DO NOT** immerse the unit for cleaning.

DO NOT operate the unit without a vessel/sample on the top plate.

DO NOT operate the unit if it shows signs of electrical or mechanical damage.

Earth Ground - Protective Conductor Terminal

Alternating Current

STANDARDS & REGULATIONS

VWR international hereby declares under its sole responsibility that the products comply with the requirements of the following directives and associated standards:

Associated EU directives:

EMC directive	2004/108/EC
LVD directive	2006/95/EC
ROHS directive	2002/95/EC
WEEE directive	2002/96/EC

Safety standards:

EN 61010-1
EN 61010-2-010
UL 61010-1

EMC standards:

EN 45501	EN 61000-4-3
EN 55022A	EN 61000-4-4
EN 61326	EN 61000-4-5
EN 6100-3-2	EN 61000-4-6
EN 61000-3-3	EN 61000-4-11
EN 61000-4-2	

SPECIFICATIONS

Overall dimensions (L x W x H):	16,5 x 14,6 x 12,1cm
Top plate dimensions (dia.):	11,4cm
Top plate material:	aluminum
Electrical (50/60 Hz):	230 volts, 0,25 amps, 20 watts
Fuses:	5mm x 20mm, 5 amp quick acting
Speed range:	100 to 1200rpm
Speed stability:	+/-2%
Weight capacity:	1000mL
Controls:	stir knob, 1 to 10 dial markings
Ship weight:	1,6kg
European Catalogue Numbers:	Euro plug: 444-2856
	UK plug: 444-2857
	Swiss plug: 444-2858

OPERATING INSTRUCTIONS

The Mini Auto Stirrer has a hinged support plate that provides automatic on-off control, thus allowing the speed control setting to remain undisturbed during each batch operation. The plate activates a motor controlled micro-switch when under the weight of a stirring vessel. The activation weight is approximately 50 grams.

1. **Getting ready:** The stir knob has a built in on-off switch at its extreme counter-clockwise position. Turn the stir knob to the off position. Plug the power cord into a properly grounded outlet. Place a vessel with solution and the appropriate accessories in the center of the top plate. This is important for stirring because the stir bar will be well within the magnetic field.
2. **Turning unit on:** From the off position, turn the stir knob clockwise until a click is heard, then continue turning until the desired speed is obtained. The stirrer may initially have to be set at a higher speed to overcome friction in the motor and the material being stirred. Once this is done, the stirrer will maintain the approximate speed set for as long as desired.
3. **Setting speed:** Speed adjustments can be made without interrupting stirring by turning the stir knob clockwise to increase speed or counter-clockwise to decrease speed.
4. **Stop stirring:** When the stirring is complete, lift the vessel off the hinged top plate. The motor will shut off. The unit is ready to stir the next sample. Change the speed setting only when needed.

OPERATING TIPS

The stirrer increases speed at a steady rate until the set-point is reached, if the stir bar is too large or the liquid is too viscous, the stirrer may not reach its set-point. The set-point speed needs to be reduced.

Stirring vessels in oil baths: When stirring a reaction vessel within an oil bath or similar set-up, the stirring function will stir up to approximately 2,54cm from the top plate. The stirring speed will vary according to liquid viscosity, stir bar length and distance from the top plate. Adjust one or all of these to achieve the desired stirring speed.

EXAMPLE: The closer the reaction vessel is to the top plate the stronger the magnetic connection.

TECHNICAL SERVICE

For more information or technical assistance contact your local VWR representative or visit www.vwr.com.

Web resources: Visit VWR's website at www.vwr.com for:

- Complete technical service contact information
- Access to VWR's Online Catalogue, and information about accessories and related products
- Additional product information and special offers

TROUBLESHOOTING

Problem	Cause	Solution
Unit will not reach desired speed	Sample volume or viscosity	Reduce volume or viscosity. If problem persists, please contact your VWR representative for repair.
Spin bar decouples	Faulty spin bar	Replace spin bar or use alternate type.
Unit is excessively noisy	Motor	This error cannot be fixed by the end user. Please contact your VWR representative for repair.

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hereby declare that the products, mentioned in Schedule comply with the essential requirements of the following Directives of the European Parliament and council:

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Title of Council Directive on harmonization of the laws of the member states 73/23/EEC of 19 February 1973 on the harmonisation of the laws of Member States relating to Electrical Equipment designed for use within certain voltage limits
89/336/EEC of 3 May 1989 on the approximation of the laws of the Member States relating to Electromagnetic Compatibility

and are in conformity with the following standard(s) and/or other normative document(s):

et sont conformes à la (aux) norme(s) ou autre(s) document(s) normatif(s) :
und der (die) folgende(n) Norm(en) oder ander(er) normativ(er) Dokument(e) entsprechen:

Reference number	Title of the applied standards:
EN 61010-2-010	Safety requirements for electrical equipment for measurement, control, and laboratory use — Part 2-010: Particular requirements for laboratory equipment for the heating of material
EN 61010-1	Safety requirements for electrical equipment for measurement, control, and laboratory use — Part 1: General requirements
EN 45501	Relating to Electromagnetic compatibility
EN55022A	
EN 61000-3-2,-3,-4-2,-4-3,-4-4,-4-5,-4-6,-4-11	
EN 61326	

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Yves Van Damme
European Regulatory Affairs Manager

Frankie Vangeel
Managing Director

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Ireland / Northern Ireland

VWR International Ltd / VWR International
(Northern Ireland) Ltd
Orion Business Campus
Northwest Business Park
Ballycoolin
Dublin 15
Tel.: 01 88 22 222
Fax: 01 88 22 333
E-mail: sales@ie.vwr.com

Italy

VWR International s.r.l.
Via Stephenson 94
20157 Milano (MI)
Tel.: 02 332 03 11
Fax: 800 152 999
E-mail: info@it.vwr.com

The Netherlands

VWR International B.V.
Postbus 8198
1005 AD Amsterdam
Tel.: 020 4808 400
Fax: 020 4808 480
E-mail: info@nl.vwr.com

Norway

VWR International AS
Haavard Martinsens vei 30
0978 Oslo
Tel.: 02290
Fax: 815 00 940
E-mail: info@no.vwr.com

Poland

Labart Sp. z o.o.
A VWR International Company
Limbowa 5
80-175 Gdansk
Tel.: 058 32 38 210
Fax: 058 32 38 205
E-mail: labart@labart.pl

Portugal

VWR International - Material de
Laboratório, Lda
Edifício Neopark
Av. Tjomas Ribeiro, 43-3 D
2790-221 Carnaxide
Tel.: 21 3600 770
Fax: 21 3600 798/9
E-mail: info@pt.vwr.com

Singapore

VWR Singapore Pte Ltd
18 Gul Drive
Singapore 629468
Tel.: +65 6505 0760
Fax: +65 6264 3780
E-mail: sales@sg.vwr.com

Spain

VWR International Eurolab S.L.
C/Tecnologia 5-17
A-7 Llinars Park
08450 - Llinars del Vallés
Barcelona
Tel.: 902 222 897
Fax: 902 430 657
E-mail: info@es.vwr.com

Sweden

VWR International AB
Fagerstagatan 18a
163 94 Stockholm
Tel.: 08 621 34 00
Fax: 08 621 34 66
E-mail: info@se.vwr.com

Switzerland

VWR International AG
Lerzenstrasse 16/18
8953 Dietikon
Tel.: 044 745 13 13
Fax: 044 745 13 10
E-mail: info@ch.vwr.com

UK

VWR International Ltd
Customer Service Center
Hunter Boulevard
Magna Park
Lutterworth
Leicestershire
LE17 4XN
Tel.: 0800 22 33 44
Fax: 01455 55 85 86
E-mail: uksales@uk.vwr.com

Legal Manufacturer:

VWR International bvba • Researchpark Haasrode 2020 • Geldenaaksebaan 464 • B-3001 Leuven • + 32 16 385011 • <http://be.vwr.com>
715088-00 (Rev 2)